

## **Materials Testing**



## SM1002

## Bench Top Tensile Testing Machine

A laboratory-scale hand-driven bench top tensile testing machine, 20 kN capacity





- Bench Top Tensile Testing Machine designed for student use
- Simple hand-operated load application for safe and easy operation
- · Digital displays of load and overall displacement
- Supplied with chucks for standard 20 mm<sup>2</sup> specimens
- Complements TecQuipment's Torsion Testing Machine (SM1001)
- Optional Extensometer (SM1002A) for Young's Modulus tests
- Optional Compression Cage and Brinell Test Set (SM1002B and SM1002C)
- Works with TecQuipment's Versatile Data Acquisition System (VDAS®)
- TecQuipment Ltd, Bonsall Street, Long Eaton, Nottingham NG10 2AN, UK
- T +44 115 972 2611 F +44 115 973 1520 E info@tecquipment.com W www.tecquipment.com
- An ISO 9001 certified company
- VDAS is a registered trademark of TecQuipment Ltd



### SM1002

## Bench Top Tensile Testing Machine

#### **Description**

A small-scale machine that fits on a benchtop and allows simple tensile tests of metal specimens up to a maximum load of 20 kN. This machine is a good partner to TecQuipment's Torsion Testing Machine (SM1001).

The machine has an extruded aluminium bed that holds the load application and load measuring mechanisms. 'Tie bars' add rigidity to the structure.

The load application mechanism includes a hand-driven worm-and-wheel gearbox, driving a lead screw with approximately 400 mm of travel. The mechanism uses ball races and self-aligning ball thrust races in the direction of loading. These low-friction bearings, with the large handwheel allow the user to apply maximum load with minimum effort. They also give smooth and progressive operation, necessary to help the user apply a steady strain rate for best results.

The unit also has a smaller "quick advance" handwheel that allows the user to set the distance between the chucks simply and quickly before each test.

The load measuring mechanism is a strain-gauged load cell that connects to a microprocessor-controlled digital display. The load display unit has a 'peak hold' function to register the maximum load before the specimen breaks.

A sliding digital display measures the tensile displacement (extension) over the entire movement. An optional precision extensometer (SM1002A) is available for increased strain measurement accuracy to allow measurements of the material's Young's Modulus.

Both the load, extension and extensometer displays can connect to TecQuipment's optional VDAS®.

The tensile specimens mount between the load application mechanism and load cell, in collet chucks via ball-jointed spigots. This ensures purely axial loading.

The equipment includes collet chucks to fit both the long and short style of TecQuipment's 20mm<sup>2</sup> specimens.

TecQuipment supply a starter set of tensile specimens with the machine, made of two different carbon steel alloys (each in their 'as drawn' and their annealed condition), brass and aluminium. TecQuipment can also supply extra specimens (contact our sales team for details). The User Guide includes drawings to allow the user to make their own specimens if needed.

For quick and reliable tests, TecQuipment's optional VDAS® gives accurate real-time data capture, monitoring and display, calculation and charting of all important readings on a computer (computer not included).

#### **Standard Features**

- Supplied with comprehensive user guide
- Five-year warranty
- Manufactured in accordance with the latest European Union directives

#### **Experiments**

- Tensile tests up to 20 kN on specimens made of different metals, to find material characteristics such as upper and lower yield strengths, tensile strength and overall extension.
- Tests of Young's Modulus (E) for the specimen material (needs SM1002A and TL specimens).

#### **Available Experiment Modules**

Brinell Test Set (SM1002C)

#### **Recommended Ancillaries**

- Bench-mounted version of the Versatile Data Acquisition System (VDAS-B)
- Extensometer (SM1002A)
- Extra TS and TL specimens refer to separate datasheet

#### **Essential Services**

Bench space needed: Approximately 1.2 m x 0.5 m

Electrical Supply:

Single-phase 100 VAC to 240 VAC, 50 Hz to 60 Hz with earth

#### **Operating Conditions**

*Operating environment:* Laboratory

Storage temperature range:

-25°C to +55°C (when packed for transport)

Operating temperature range:

+5°C to +40°C

Operating relative humidity range:

80% at temperatures <  $31^{\circ}$ C decreasing linearly to 50% at  $40^{\circ}$ C

#### **Sound Levels**

Less than 70 dB(A)



- T +44 115 972 2611 F +44 115 973 1520 E info@tecquipment.com W www.tecquipment.com
- An ISO 9001 certified company
- VDAS is a registered trademark of TecQuipment Ltd



## **SM1002**

## Bench Top Tensile Testing Machine

#### **Specifications (SM1002)**

Nett dimensions and weight: 1100 mm long x 350 mm high x 400 mm wide and 41.5 kg (including display unit)

Approximate packed volume and weight: 0.3 m<sup>3</sup> and 80 kg

#### Digital Displays:

- Load display resolution of 0.01 kN
- Extension display resolution of 0.01 mm

Tensile Test specimens included (30 in sets of 5):

- 0.1% and 0.4% carbon steel (as drawn)
- 0.1% and 0.4% carbon steel (normalised)
- Brass and Aluminium

Each set of five specimens has  $2 \times long$  (TL) specimens and  $3 \times long$  (TS) specimens.

#### Specifications (SM1002A)

Nett dimensions and weight: 120 mm x 100 mm x 65 mm and 1 kg

Approximate packed volume and weight: 0.005 m<sup>3</sup> and 2 kg

Display resolution of 0.001 mm



- T +44 115 972 2611 F +44 115 973 1520 E info@tecquipment.com W www.tecquipment.com
- An ISO 9001 certified company
- VDAS is a registered trademark of TecQuipment Ltd



# tradition. innovation. integration.

**infoWERK** is a leading expert in the development of eLearning courseware, learning system solutions, teaching and AV equipment.

Furthermore infoWERK is the representative and system integrator of "TecQuipment".

**TecQuipment** is one of the global leaders in technical teaching equipment for engineering. If you are interested in one of TecQuipment's products feel free to contact us at:



E-Mail: info@de.infowerk.at / Website: infowerk.at